碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301或800-8283301
订货e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

Ad-GFP-LC3B

产品编号	产品名称	包装
C3006-1ml	Ad-GFP-LC3B	1ml
C3006-10ml	Ad-GFP-LC3B	10ml

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Species	Cell line	Origin	Transduction efficiency
Human	293	Embryonic Kidney	95%
	MCF-10A	Mammary epithelium	51%
	MDA-MB-231	Malignant breast cancer cell	72%
	HepG2	Hepatocarcinoma	70%
	T24	Transitional cell carcinoma	37%
	TCCSUP	Transitional cell carcinoma	53%
	PC-3	Prostate cancer metastases site	67%
	DU145	Prostate cancer metastases site	86%
	LNCaP	Prostate cancer metastases site	69%
	Jurkat	T cell leukemia	2%
	HRT	Rectal tumor	5%
Murine	MT1A2	Mammary tumor transformed with SV40 T antigen	92%
	NIH 3T3	Mouse fibroblast	12%
	WEHI	Fibrosarcoma	10%
	AML12	Liver	28%
	MS-K68	Spleen	48%
	EL4	T cell	2%
Bovine	FBRT HE1	Fetal retina transformed with Had5 E1	90%
	MDBK	Kidney	70%
	FBK-34	Kidney transformed with BAd3 E1	25%
	BT-5705	Turbinate	28%
	EBL	Embryonic lung	8%
Porcine	FPRT HE1-5	Fetal retina transformed with HAd5 E1	91%
	PK-15	Kidney	76%
	PK-M12	Kidney cell transformed with SV40 T antigen	71%
	PT-K75	Turbinate	22%
	EPL	Embryonic lung	53%
	SBC-1765	Derived from buffy coat cells	73%

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Table 1. Summary of Recommended Biosafety Levels for Infectious Agents

BSL	Agents	Practices	Primary Barriers and Safety Equipment	Facilities (Secondary Barriers)
1	Not known to consistently cause diseases in healthy adults	Standard microbiological practices	<ul style="list-style-type: none"> ■ No primary barriers required. ■ PPE: laboratory coats and gloves; eye, face protection, as needed 	Laboratory bench and sink required
2	<ul style="list-style-type: none"> ■ Agents associated with human disease ■ Routes of transmission include percutaneous injury, ingestion, mucous membrane exposure 	BSL-1 practice plus: <ul style="list-style-type: none"> ■ Limited access ■ Biohazard warning signs ■ “Sharps” precautions ■ Biosafety manual defining any needed waste decontamination 	Primary barriers: <ul style="list-style-type: none"> ■ BSCs or other physical containment devices used for all manipulations of agents that cause splashes or aerosols of infectious materials ■ PPE: Laboratory coats, 	BSL-1 plus: <ul style="list-style-type: none"> ■ Autoclave available

		or medical surveillance policies	gloves, face and eye protection, as needed	
3	Indigenous or exotic agents that may cause serious or potentially lethal disease through the inhalation route of exposure	BSL-2 practice plus: <ul style="list-style-type: none"> ■ Controlled access ■ Decontamination of all waste ■ Decontamination of laboratory clothing before laundering 	Primary barriers: <ul style="list-style-type: none"> ■ BSCs or other physical containment devices used for all open manipulations of agents ■ PPE: Protective laboratory clothing, gloves, face, eye and respiratory protection, as needed 	BSL-2 plus: <ul style="list-style-type: none"> ■ Physical separation from access corridors ■ Self-closing, double-door access ■ Exhausted air not recirculated ■ Negative airflow into laboratory ■ Entry through airlock or anteroom ■ Hand washing sink near laboratory exit
4	<ul style="list-style-type: none"> ■ Dangerous/exotic agents which pose high individual risk of aerosol-transmitted laboratory infections that are frequently fatal, for which there are no vaccines or treatments ■ Agents with a close or identical antigenic relationship to an agent requiring BSL-4 until data are available to redesignate the level ■ Related agents with unknown risk of transmission 	BSL-3 practices plus: <ul style="list-style-type: none"> ■ Clothing change before entering ■ Shower on exit ■ All material decontaminated on exit from facility 	Primary barriers: <ul style="list-style-type: none"> ■ All procedures conducted in Class III BSCs or Class I or II BSCs in combination with full-body, air-supplied, positive pressure suit 	BSL-3 plus: <ul style="list-style-type: none"> ■ Separate building or isolated zone ■ Dedicated supply and exhaust, vacuum, and decontamination systems ■ Other requirements outlined in the text

BSL, biosafety level; PPE, personal protective equipment.

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产品编号	产品名称	包装
C3006-1ml	Ad-GFP-LC3B	1ml
C3006-10ml	Ad-GFP-LC3B	10ml
C3011-1ml	Ad-mCherry-GFP-LC3B	1ml
C3010-10ml	Ad-mCherry-GFP-LC3B	10ml
C3012-1ml	AdPlus-mCherry-GFP-LC3B	1ml
C3012-10ml	AdPlus-mCherry-GFP-LC3B	10ml

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- Zhao W, Peng F, Shu M, Liu H, Hou X, Wang X, Ye J, Zhao B, Wang K, Zhong C, Xue L, Gao M, Liu Y, Zhao S. Isogambogenic Acid Inhibits the Growth of Glioma Through Activation of the AMPK-mTOR Pathway. *CELL PHYSIOL BIOCHEM*. 2017;44(4):1381-1395.
- Zhao Q, Niu Q, Chen J, Xia T, Zhou G, Li P, Dong L, Xu C, Tian Z, Luo C, Liu L, Zhang S, Wang A. Roles of mitochondrial fission inhibition in developmental fluoride neurotoxicity: mechanisms of action in vitro and associations with cognition in rats and children. *Arch Toxicol*. 2019 Mar;93(3):709-726.
- Fu XT, Song K, Zhou J, Shi YH, Liu WR, Shi GM, Gao Q, Wang XY, Ding ZB, Fan J. Tumor-associated macrophages modulate resistance to oxaliplatin via inducing autophagy in hepatocellular carcinoma. *Cancer Cell Int*. 2019 Mar 25;19:71.
- Chu P, He L, Yang C, Li Y, Huang R, Liao L, Li Y, Zhu Z, Wang Y. Characterisation and function of TRIM23 in grass carp (*Ctenopharyngodon idella*). *Fish Shellfish Immunol*. 2019 May;88:627-635.
- Chu P, He L, Yang C, Zeng W, Huang R, Liao L, Li Y, Zhu Z, Wang Y. Grass carp ATG5 and ATG12 promote autophagy but down-regulate the transcriptional expression levels of IFN- γ signaling pathway. *Fish Shellfish Immunol*. 2019 Sep;92:600-611.
- Shi X, Zhang H, Jin W, Liu W, Yin H, Li Y, Dong H. Metronomic photodynamic therapy with 5-aminolevulinic acid induces apoptosis and autophagy in human SW837 colorectal cancer cells. *J Photochem Photobiol B*. 2019 Sep;198:111586.
- Wang P, Zhang SD, Jiao J, Wang W, Yu L, Zhao XL, Wang LT, Meng D, Fu YJ. ROS-mediated p53 activation by juglone enhances apoptosis and autophagy in vivo and in vitro. *Toxicol Appl Pharmacol*. 2019 Sep 15;379:114647.
- Chu P, He L, Yang C, Zeng W, Huang R, Liao L, Li Y, Zhu Z, Wang Y. Grass carp ATG5 and ATG12 promote autophagy but down-regulate the transcriptional expression levels of IFN- γ signaling pathway. *Fish Shellfish Immunol*. 2019 Sep;92:600-611.
- Shi X, Zhang H, Jin W, Liu W, Yin H, Li Y, Dong H. Metronomic photodynamic therapy with 5-aminolevulinic acid induces apoptosis and autophagy in human SW837 colorectal cancer cells. *J Photochem Photobiol B*. 2019 Sep;198:111586.
- Wang P, Zhang SD, Jiao J, Wang W, Yu L, Zhao XL, Wang LT, Meng D, Fu YJ. ROS-mediated p53 activation by juglone enhances apoptosis and autophagy in vivo and in vitro. *Toxicol Appl Pharmacol*. 2019 Sep 15;379:114647.

11. Tan X,Zou L,Qin J,Xia D,Zhou Y,Jin G,Jiang Z,Li H.SQSTM1/p62 is involved in docosahexaenoic acid-induced cellular autophagy in glioblastoma cell lines.In *Vitro Cell Dev Biol Anim.* 2019 Oct;55(9):703-712.
12. Tang W,Lin D,Chen M,Li Z,Zhang W,Hu W,Li F.PTEN-mediated mitophagy and APE1 overexpression protects against cardiac hypoxia/reoxygenation injury.In *Vitro Cell Dev Biol Anim.* 2019 Oct;55(9):741-748.
13. Huang Y,Li Y,Qu Y,Zheng Y,Ouyang M,Zhang Y,Lai W,Xu Q.UVA-induced photoaging inhibits autophagic degradation by impairing lysosomal function in dermal fibroblasts.*Biochem Biophys Res Commun.* 2019 Oct 22;518(4):611-618.
14. Danyao Nie,Xinhua Liu,Yuan Wang,Wenling He,Ming Li,Yun Peng,Jing Zhang,Liangnan Sun,Zonghui Yan,Lin Ye.Involvement of the JNK signaling in granular corneal dystrophy by modulating TGF- β -induced TGFBI expression and corneal fibroblast apoptosis IN *VITRO CELL DEV-AN.* 2020 Mar;56(3):234-242.;doi: 10.1007/s11626-019-00424-6
15. Na Geng,Xiaozhou Wang,Xiaohui Yu,Run Wang,Yiran Zhu,Meihua Zhang,Jianzhu Liu,Yongxia Liu.Staphylococcus aureus Avoids Autophagy Clearance of Bovine Mammary Epithelial Cells by Impairing Lysosomal Function *Front Immunol.* 2020 May 5;11:746.;doi: 10.3389/fimmu.2020.00746
16. Yun Tsai,Changbo Xia,Zhongwen Sun.The Inhibitory Effect of 6-Gingerol on Ubiquitin-Specific Peptidase 14 Enhances Autophagy-Dependent Ferroptosis and Anti-Tumor in vivo and in vitro *Front Pharmacol.* 2020 Nov 13;11:598555.;doi: 10.3389/fphar.2020.598555
17. Ping Zhang,Hongfei Ni,Yang Zhang,Wenping Xu,Jufang Gao,JIAGAO Cheng,Liming Tao.Ivermectin confers its cytotoxic effects by inducing AMPK/mTOR-mediated autophagy and DNA damage *Chemosphere.* 2020 Nov;259:127448.;doi: 10.1016/j.chemosphere.2020.127448
18. Shuai Wang,Rong Guo,Yongxia Su,Chen Yang,Yazhou Guo,Chengjian Tan,Baoyu Zhao.Swainsonine promotes apoptosis by impairing lysosomal function and inhibiting autophagic degradation in rat primary renal tubular epithelial cells *CHEM-BIOL INTERACT.* 2021 Feb 25;336:109319.;doi: 10.1016/j.cbi.2020.109319

Version 2021.09.01