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1. Jia Y,Liu X.Expression of Na⁺/K⁺-ATPase Was Affected by Salinity Change in Pacific abalone *Haliotis discus hannai*.*Front Physiol* . 2018 Sep 7;9:1244.
2. Gu Z,Zhao L,Ge C,Liu S,Fang G,Chen SS,Yang Z,Zhou R.Facet-regulated adhesion of double-stranded DNA on palladium surfaces.*Nanoscale*. 2019 Jan 23;11(4):1827-1836.
3. Yang L,Sun J,Liu Q,Zhu R,Yang Q,Hua J,Zheng L,Li K,Wang S,Li A.Synergetic Functional Nanocomposites Enhance Immunotherapy in Solid Tumors by Remodeling the Immunoenvironment.*Adv Sci (Weinh)*. 2019 Feb 16;6(8):1802012.
4. Ji X,Tang Z,Shuai W,Zhang Z,Li J,Chen L,Cao J,Yin W.Endogenous peptide LYENRL prevents the activation of hypertrophic scar-derived fibroblasts by inhibiting the TGF-β1/Smad pathway.*Life Sci*. 2019 Aug 15;231:116674.
5. Peng T,Deng X,Tian F,Li Z,Jiang P,Zhao X,Chen G,Chen Y,Zheng P,Li D,Wang S.The interaction of LOXL2 with GATA6 induces VEGFA expression and angiogenesis in cholangiocarcinoma.*Int J Oncol*. 2019 Sep;55(3):657-670.
6. Hang Gao,Siyue Lou,Huanwu Hong,Qiufu Ge,Huajun Zhao.Autophagy Inhibition Potentiates the Anticancer Effects of a Bendamustine Derivative NL-101 in Acute T Lymphocytic Leukemia *Biomed Res Int*. 2020 Feb 14;2020:1520651.;doi: 10.1155/2020/1520651
7. Mengzhu Wang,Zhigang Wang,Bin Qiao,Jin Cao,Luya Quan,Yuanli Luo,Hanwen Qi,Xiaowen Zhong,Yubei He,Xianquan Zhang,Lan Hao.Inhibited Metastasis and Amplified Chemotherapeutic Effects by Epigene-Transfection Based on a Tumor-Targeting Nanoparticle *INT J NANOMED*. 2020 Jun 23;15:4483-4500.;doi: 10.2147/IJN.S247567

Version 2021.09.01