碧云天生物技术/Beyotime Biotechnology  
订货热线: 400-168-3301 或 800-8283301  
订货 e-mail: order@beyotime.com  
技术咨询: info@beyotime.com  
网址: http://www.beyotime.com

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1. Wang C, Xie H, Liu X, Qin Q, Wu X, Liu H, Liu C. Role of nuclear factor- $\kappa$ B in volatile anaesthetic preconditioning with sevoflurane during myocardial ischaemia/reperfusion. *Eur J Anaesthesiol.* 2010 Aug;27(8):747-56.
2. Wu Y, Feng W, Zhang H, Li S, Wang D, Pan X, Hu S. Ca<sup>2+</sup>-regulatory proteins in cardiomyocytes from the right ventricle in children with congenital heart disease. *J Transl Med.* 2012 Apr 2;10:67.
3. Wang X, Lin H, Sui J, Cao L. The effect of fish matrix on the enzyme-linked immunosorbent assay of antibiotics. *J Sci Food Agric.* 2013 May;93(7):1603-9.
4. Wang HJ, Tan YZ. Methods for assessing effects of Wnt/ $\beta$ -catenin signaling in senescence of mesenchymal stem cells. *Methods Mol Biol.* 2013;976:111-30.
5. Xie H, Liu X, Wang C, Zhu J, Yang C, Liu C, Liu H, Wu X. The changes of technetium-99m-labeled annexin-V in delayed anesthetic preconditioning during myocardial ischemia/reperfusion. *Mol Biol Rep.* 2014 Jan;41(1):131-7.
6. Jiang X, Ge H, Zhou C, Chai X, Deng H. The role of transforming growth factor  $\beta$ 1 in fractional laser resurfacing with a carbon dioxide laser. *Lasers Med Sci.* 2014 Mar;29(2):681-7.
7. Sun L, Li G, Dai B, Tan W, Zhao H, Li X, Wang A. Silence of MACC1 expression by RNA interference inhibits proliferation, invasion and metastasis, and promotes apoptosis in U251 human malignant glioma cells. *Mol Med Rep.* 2015 Sep;12(3):3423-31.
8. Cui M, Lin H, Wang X, Cao L, Sui J. 5-Sulfosalicylic acid dihydrate-based pretreatment for the modification of enzyme-linked immunoassay of fluoroquinolones in fishery products. *J Immunoassay Immunochem.* 2015;36(5):517-31.
9. Jiang F, Zhang Z, Zhang Y, Wu J, Yu L, Liu S. L carnitine ameliorates the liver inflammatory response by regulating carnitine palmitoyltransferase I dependent PPAR $\gamma$  signaling. *Mol Med Rep.* 2016 Feb;13(2):1320-8.
10. Li Q, Tang Y, Qin J, Yi L, Yang Y, Wang J, He J, Su Q, Tan H. Subcellular localization of DJ-1 in human HL-60 leukemia cells in response to diallyl disulfide treatment. *Mol Med Rep.* 2016 Nov;14(5):4666-4672.
11. Lin JP, Chen CQ, Huang LE, Li NN, Yang Y, Zhu SM, Yao YX. Dexmedetomidine Attenuates Neuropathic Pain by Inhibiting P2X7R Expression and ERK Phosphorylation in Rats. *Exp Neurobiol.* 2018 Aug;27(4):267-276.
12. Wang X, Yi K, Zhao Y. Fucoidan inhibits amyloid- $\beta$ -induced toxicity in transgenic *Caenorhabditis elegans* by reducing the accumulation of amyloid- $\beta$  and decreasing the production of reactive oxygen species. *Food Funct.* 2018 Jan 24;9(1):552-560.
13. Pan R, Jiang Y, Sun L, Wang R, Zhuang K, Zhao Y, Wang H, Ali MA, Xu H, Man C. Gold nanoparticle-based enhanced lateral flow immunoassay for detection of

Cronobacter sakazakii in powdered infant formula. J Dairy Sci . 2018 May;101(5):3835-3843.

14. Liu K,He K,Xue T,Liu P,Xu LXThe cryo-thermal therapy-induced IL-6-rich acute pro-inflammatory response promoted DCs phenotypic maturation as the prerequisite to CD4+ T cell differentiation. INT J HYPERTHER . 2018 May;34(3):261-272.
15. Wang T,Hu J,Luo H,Li H,Zhou J,Zhou L,Wei SPhotosensitizer and Autophagy Promoter Coloaded ROS-Responsive Dendrimer-Assembled Carrier for Synergistic Enhancement of Tumor Growth Suppression. Small . 2018 Sep;14(38):e1802337.

Version 2019.10.17