碧云天生物技术/Beyotime Biotechnology
 订货热线: 400-1683301或800-8283301
 订货e-mail: order@beyotime.com
 技术咨询: info@beyotime.com
 网址: http://www.beyotime.com

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- Xin Jiang, Wenjing Chen, Hang Su, Fuguo Shen, Wenlong Xiao, Wencai Sun Puerarin facilitates osteogenesis in steroid-induced necrosis of rabbit femoral head and osteogenesis of steroid-induced osteocytes via miR-34a upregulation Cytokine 2021 Jul;143:155512.
- Hong Wang, Shuqin Wang, Shanshan Huang MiR-494-3p alleviates acute lung injury through regulating NLRP3 activation by targeting CMPK2 Biochem Cell Biol 2021 Jun;99(3):286-295.
- Weilan Wang, Chengmin Huang, Peng Luo, Jiang Yao, Jie Li, Wenxia Wang, Fengqin Liu Circular RNA circWDR27 Promotes Papillary Thyroid Cancer Progression by Regulating miR-215-5p/TRIM44 Axis Onco Targets Ther 2021 May 19;14:3281-3293.
- Xiong Qian, Qi Xiao, Zongqi Li Tectorigenin regulates migration, invasion, and apoptosis in dexamethasone-induced human airway epithelial cells through up-regulating miR-222-3p Drug Dev Res 2021 Nov;82(7):959-968.
- Chen Shao, Guoqi Niu, Peng Su, Jingquan Zhang, Xunbing Zhu, Guansheng Han, Panpan Xu, Jianzhong Bai, Kui Sun, Yongming Sun CircFOXK2 Promotes the Progression of Osteoarthritis by Regulating the miR-4640-5p/NOTCH2 Axis Mod Rheumatol 2022 Dec 20:roac158.
- Bo Feng, Ke Chen, Weiwei Zhang, Qi Zheng, Yong He Silencing of lncRNA MIR31HG promotes nasopharyngeal carcinoma cell proliferation and inhibits apoptosis through suppressing the PI3K/AKT signaling pathway J Clin Lab Anal 2022 Dec;36(12):e24720.
- Hang Li, Lixin Chai, Zujun Ding, Huabo He CircCOL1A2 Sponges MiR-1286 to Promote Cell Invasion and Migration of Gastric Cancer by Elevating Expression of USP10 to Downregulate RFC2 Ubiquitination Level J Microbiol Biotechnol 2022 Jul 28;32(7):938-948.
- Liming Cao, Fan Lin TECPR1 Induces Apoptosis in Non-Small Cell Lung Carcinoma via ATG5 Upregulation-Induced Autophagy Promotion Ann Clin Lab Sci 2022 Jul;52(4):580-592.
- Wei Li, Kai Li, Zhong Wang, Zhenzhong Fa MicroRNA-377-3p promotes cell proliferation and inhibits cell cycle arrest and cell apoptosis in hepatocellular carcinoma by affecting EGR1-mediated p53 activation Pathol Res Pract 2022 Jun;234:153855.
- Hong Lu, Li Ai, Baoyue Zhang TNF-α induces AQP4 overexpression in astrocytes through the NF-κB pathway causing cellular edema and apoptosis Biosci Rep 2022 Mar 31;42(3):BSR20212224.
- Weitao Yao, Jingyu Hou, Guoqing Liu, Fangxing Wu, Qiang Yan, Liangyu Guo, Chuchu Wang LncRNA CBR3-AS1 promotes osteosarcoma progression through the network of miR-140-5p/DDX54-NUCKS1-mTOR signaling pathway Mol Ther Oncolytics 2022 Mar 8;25:189-200.

12. Cuicui Wu β -catenin inhibitors ICG-001 and pyrvinium sensitize bortezomib-resistant multiple myeloma cells to bortezomib *Oncol Lett* 2022 May 12;24(1):205.
13. Yanbo Zhu, Fang Li, Yilong Wan, Hansi Liang, Si Li, Bo Peng, Liqun Shao, Yunyun Xu, Dong Jiang Cancer-Secreted Exosomal MiR-620 Inhibits ESCC Aerobic Glycolysis via FOXM1/HER2 Pathway and Promotes Metastasis *Front Oncol* 2022 May 16;12:756109.
14. Yuli Ge, Huiling Sun, Lanman Xu, Weiping Zhang, Jiaojian Lv, Yongping Chen The amelioration of alcohol-induced liver and intestinal barrier injury by *Lactobacillus rhamnosus* Gorbach-Goldin (LGG) is dependent on Interleukin 22 (IL-22) expression *Bioengineered* 2022 May;13(5):12650-12660.
15. Jing Yuan, Lili Wang, Shuang Han, Zhongli Wang, Yanhui Ni, Yanping Geng, Li Zhang PPAR γ Silencing Improves Blood Pressure Control and Alleviates Renal Damage by Modulating RAS Circadian Rhythm in Hypertensive Rats *Ann Clin Lab Sci* 2022 May;52(3):452-461.
16. Guangzhi Li, Dongmei Zhang, Shizhen Wang, Ni Jie, Yongting Qin Protective Role of Taraxasterol against Cardiovascular Aging and Aging-Induced Desensitization of Insulin Signaling *Front Biosci (Landmark Ed)* 2022 Nov 22;27(11):311.
17. Yuehong Wang, Zhilian Wang, Keyan Cheng, Qirong Hao FAM201A Promotes Cervical Cancer Progression and Metastasis through miR-1271-5p/Flotillin-1 Axis Targeting-Induced Wnt/ β -Catenin Pathway *J Oncol* 2022 Oct 3:2022:1123839.
18. Yi Liu, Shanshan Ma, Qichao Ma, Haigang Zhu Silencing LINC00665 inhibits cutaneous melanoma in vitro progression and induces apoptosis via the miR-339-3p/TUBB *J Clin Lab Anal* 2022 Sep;36(9):e24630.
19. Yong Zhang, Yuan Fang, Lijie Ma, Jing Xu, Chentao Lv, Li Deng, Guanghui Zhu LINC00857 regulated by ZNF460 enhances the expression of CLDN12 by sponging miR-150-5p and recruiting SRSF1 for alternative splicing to promote epithelial-mesenchymal transformation of pancreatic aden *RNA Biol* 2022;19(1):548-559.
20. Lantao Zhao, Hongya Xie, Peng Li, Haisheng Chen, Jiang He, Liang Wang, Yonghong Wang, Bin Ni CircTFF1 Promotes Proliferation, Migration and Invasion of Lung Cancer Cells by Facilitating Methylation of BCL6B Promoter via miR-29c-3p/DNMT3A Axis *Mol Biotechnol* 2023 Jun;65(6):942-952.

Version 2024.03.12