碧云天生物技术/Beyotime Biotechnology
 订货热线: 400-1683301或800-8283301
 订货e-mail: order@beyotime.com
 技术咨询: info@beyotime.com
 网址: http://www.beyotime.com

Anti-Myc Affinity Gel (Anti-Myc亲和凝胶)

| 产品编号 | 产品名称 | 包装 |
|-------------|--------------------------------------|-------|
| P2285-0.5ml | Anti-Myc Affinity Gel (Anti-Myc亲和凝胶) | 0.5ml |
| P2285-2ml | Anti-Myc Affinity Gel (Anti-Myc亲和凝胶) | 2ml |
| P2285-10ml | Anti-Myc Affinity Gel (Anti-Myc亲和凝胶) | 10ml |

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| | |
|------------------------|---|
| Cat. No. | P2285 |
| Product name | Anti-Myc Affinity Gel (Anti-Myc亲和凝胶) |
| Product content | 50% settled gel in 50% glycerol with 10mM phosphate-buffered saline and preservative (pH7.4) |
| Matrix | 4% agarose |
| Average bead size | ~90µm |
| Antibody | Mouse monoclonal antibody against Myc-tag |
| Isotype | IgG1 |
| M.W. of antibody | Approximately 150kDa |
| Antibody concentration | Approximately 8mg Myc antibody per ml settled gel |
| Binding capacity | Approximately 1mg Myc-tagged protein per ml settled gel |
| Elution method | Acid, alkaline, neutral, peptide competitive or SDS-PAGE loading buffer elution. Note: If elute with SDS-PAGE loading buffer, the light (~25kDa) and heavy (~50kDa) chain of antibody will be denatured and release from the gel. |
| Reagent compatibility | Chaotropic reagents will denature the target Myc-tagged protein. Do not exceed 0.3M GuHCl or 1.5M Urea. |
| Application | Suitable for IP、Co-IP and protein purification. |
| Storage | -20°C |

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| Problem | Possible Causes | Solution |
|---|---|---|
| Large amount of tagged protein found in the flow through. | Binding time is not enough. | If using batch method, increase the binding time experimentally; If using column method, use a lower flow rate when loading samples. |
| | Column is overloaded. | Reduce the amount of the sample added to the gel or increase the amount of gel. |
| | Tag is not accessible to gel. | Expose the epitope tag by adding low amount of denaturant to the protein extract (dialysis may be needed before applying onto gel), or fuse the tag to the other terminus of the target protein. |
| | Gel has not been regenerated since last purification. | Perform gel regeneration procedure prior to binding. |
| | Reagent compatibility problem. | Dialyze the sample against TBS before purification procedure. |
| | The target protein has been degraded. | 1. Prepare fresh lysates. Avoid using frozen lysates. 2. Perform purification at lower temperature, such as 4°C. 3. Use appropriate protease inhibitors in the lysate or increase their concentrations to prevent degradation of the fusion protein. |
| Very few or no tagged protein exists in the eluate. | Protein is not completely eluted. | Change elution methods. |
| | No target protein expressed. | Make sure the protein of interest contains the tag by Western blot or dot blot analyses. |
| | Very low protein expression level. | 1. Use larger volume of cell lysate. 2. Optimize expression conditions to raise the protein expression level. |
| | Washes are too stringent. | 1. Reduce the number of washes. 2. Avoid adding high concentrations of NaCl to the mixture. 3. Use solutions that contain less or no detergent |
| | Incubation times are inadequate. | Increase the incubation times with the affinity gel (from several hours to overnight). |
| | Interfering substance is present in sample. | 1. Lysates containing high concentrations of DTT, 2-mercaptoethanol, or other reducing agents may destroy antibody function, and must be avoided. 2. Excessive detergent concentrations may interfere with the antibody-antigen interaction. Detergent levels in buffers may be reduced by dilution. |
| | Detection system is inadequate. | If Western blot detection is used: 1. Check primary and secondary antibodies using proper controls to confirm binding and reactivity. 2. Verify that the transfer was adequate by staining the membrane with Ponceau S. 3. Use fresh detection substrate or try a different detection system. |
| Multiple protein bands found in | The protein is not stable at room temperature. | Purify the target protein at lower temperature, such as 4°C. |

| | | |
|-------------------------|--|---|
| the eluate. | Protein degradation due to proteases activity during purification process. | Add protease inhibitors to cell lysate. |
| | Non-specific binding. | 1. Prepare cell lysate again. 2. Add additional wash steps. |
| Background is too high. | Proteins bind nonspecifically to the monoclonal antibody, the gel beads, or the microcentrifuge tubes. | 1. Pre-clear lysate with Mouse IgG-Agarose (P2055+A7028) to remove nonspecific binding proteins. 2. After suspending beads for the final wash, transfer entire sample to a clean microcentrifuge tube before centrifugation. |
| | Washes are insufficient. | 1. Increase the number of washes. 2. Prolong duration of the washes, incubating each wash for at least 15 minutes. 3. Increase the salt and/or detergent concentrations in the wash solutions. 4. Centrifuge at lower speed to avoid nonspecific trapping of denatured proteins from the lysate during the initial centrifugation of the affinity gel complexes. |

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| 产品编号 | 产品名称 | 包装 |
|-------------|---|---------|
| P2006 | Protein A Agarose (Fast Flow, 进口分装) | 2ml |
| P2009 | Protein G Agarose (Fast Flow, 进口分装) | 2ml |
| P2012 | Protein A+G Agarose (Fast Flow, 进口分装) | 2ml |
| P2051-2ml | Protein A Agarose (Fast Flow, for IP) | 2ml |
| P2051-10ml | Protein A Agarose (Fast Flow, for IP) | 10ml |
| P2051-50ml | Protein A Agarose (Fast Flow, for IP) | 50ml |
| P2053-2ml | Protein G Agarose (Fast Flow, for IP) | 2ml |
| P2053-10ml | Protein G Agarose (Fast Flow, for IP) | 10ml |
| P2053-50ml | Protein G Agarose (Fast Flow, for IP) | 50ml |
| P2055-2ml | Protein A+G Agarose (Fast Flow, for IP) | 2ml |
| P2055-10ml | Protein A+G Agarose (Fast Flow, for IP) | 10ml |
| P2055-50ml | Protein A+G Agarose (Fast Flow, for IP) | 50ml |
| P2265 | Mouse IgG Agarose (小鼠IgG琼脂糖凝胶) | 1ml/5ml |
| P2267 | Rabbit IgG Agarose (兔IgG琼脂糖凝胶) | 1ml/5ml |
| P2271-0.5ml | Anti-Flag Affinity Gel (Anti-Flag亲和凝胶) | 0.5ml |
| P2271-2ml | Anti-Flag Affinity Gel (Anti-Flag亲和凝胶) | 2ml |
| P2271-10ml | Anti-Flag Affinity Gel (Anti-Flag亲和凝胶) | 10ml |
| P2282-0.5ml | Anti-Flag Affinity Gel (Anti-Flag亲和凝胶) | 0.5ml |
| P2282-2ml | Anti-Flag Affinity Gel (Anti-Flag亲和凝胶) | 2ml |
| P2282-10ml | Anti-Flag Affinity Gel (Anti-Flag亲和凝胶) | 10ml |
| P2285-0.5ml | Anti-Myc Affinity Gel (Anti-Myc亲和凝胶) | 0.5ml |
| P2285-2ml | Anti-Myc Affinity Gel (Anti-Myc亲和凝胶) | 2ml |
| P2285-10ml | Anti-Myc Affinity Gel (Anti-Myc亲和凝胶) | 10ml |
| P2287-0.5ml | Anti-HA Affinity Gel (Anti-HA亲和凝胶) | 0.5ml |
| P2287-2ml | Anti-HA Affinity Gel (Anti-HA亲和凝胶) | 2ml |
| P2287-10ml | Anti-HA Affinity Gel (Anti-HA亲和凝胶) | 10ml |
| P2289-0.5ml | Anti-V5 Affinity Gel (Anti-V5亲和凝胶) | 0.5ml |
| P2289-2ml | Anti-V5 Affinity Gel (Anti-V5亲和凝胶) | 2ml |
| P2289-10ml | Anti-V5 Affinity Gel (Anti-V5亲和凝胶) | 10ml |

Version 2021.08.18