碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301 或 800-8283301
订货 e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: http://www.beyotime.com

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1. Wang G, Liang J, Gao LR, Si ZP, Zhang XT, Liang G, Yan Y, Li K, Cheng X, Bao Y, Chuai M, Chen LG, Lu DX, Yang X. . Baicalin administration attenuates hyperglycemia-induced malformation of cardiovascular system. Cell Death Dis. 2018 Feb 14;9(2):234.
2. Wang G, Liang J, Gao LR, Si ZP, Zhang XT, Liang G, Yan Y, Li K, Cheng X, Bao Y, Chuai M, Chen LG, Lu DX, Yang X. . Baicalin administration attenuates hyperglycemia-induced malformation of cardiovascular system. Cell Death Dis. 2018 Feb 14;9(2):234.
3. Tian XQ, Ma KT, Wang XW, Wang Y, Guo ZK, Si JQ. . Effects of the Calcium-Activated Chloride Channel Inhibitors T16Ainh-A01 and CaCCinh-A01 on Cardiac Fibroblast Function. CELL PHYSIOL BIOCHEM. 2018;49(2):706-716.
4. Luo L, Sun YJ, Yang L, Huang S, Wu YJ. . Avermectin induces P-glycoprotein expression in S2 cells via the calcium/calmodulin/NF- κ B pathway. CHEM-BIOL INTERACT. 2013 Apr 25;203(2):430-9.
5. Liu LI, Cai S, Qiu G, Lin J. . Fluid shear stress enhances the cell volume decrease of osteoblast cells by increasing the expression of the CLC-3 chloride channel. Biomed Rep.. 2016 Apr;4(4):408-412.
6. Qiu LY, Duan GL, Yan YF, Li YY, Wang H, Xiao L, Liao ZP, Chen HP. . Sasanquasaponin induces increase of Cl⁻/HCO₃⁻ exchange of anion exchanger 3 and promotes intracellular Cl⁻ efflux in hypoxia/reoxygenation cardiomyocytes. Mol Med Rep. 2017 Sep;16(3):2953-2961.
7. Zhu J, Xu Y, Ren G, Hu X, Wang C, Yang Z, Li Z, Mao W, Lu D. . Tanshinone IIA Sodium sulfonate regulates antioxidant system, inflammation, and endothelial dysfunction in atherosclerosis by downregulation of CLIC1. Eur J Pharmacol. 2017 Nov 15;815:427-436.
8. Zhou C, Tang X, Xu J, Wang J, Yang Y, Chen Y, Chen L, Wang L, Zhu L, Yang H. . Opening of the CLC-3 chloride channel induced by dihydroartemisinin contributed to early apoptotic events in human poorly differentiated nasopharyngeal carcinoma cells. J Cell Biochem. 2018 Nov;119(11):9560-9572.
9. Su J, Xu T, Jiang G, Hou M, Liang M, Cheng H, Li Q. . Gambogic acid triggers apoptosis in human nasopharyngeal carcinoma CNE-2Z cells by activating volume-sensitive outwardly rectifying chloride channel. Fitoterapia. 2019 Mar;133:150-158.
10. Chen Li, LinLin Chen, YuanYuan Wang, TingTing Wang, Dong Di, Hao Zhang, HuanHuan Zhao, Xu Shen, Jun Guo. . Protein Nanoparticle-Related Osmotic Pressure Modifies Nonselective Permeability of the Blood-Brain Barrier by Increasing Membrane Fluidity Int J Nanomedicine. 2021 Mar

1;16:1663-1680.

11. Zhiqin Deng, Zicong Lin, Qing Zhong, Minqiang Lu, Huankun Fang, Jianquan Liu, Li Duan, Lixin Chen, Liwei Wang, Daping Wang, Wencui Li . Interleukin 1 beta-induced chloride currents are important in osteoarthritis onset: an in vitro study *Acta Biochim Biophys Sin (Shanghai)*. 2021 Mar 26;53(4):400-409.
12. Xiang Dai, Jun Li, XiJiang Hu, Jian Ye, WenQian Cai . SLC26A4 Mutation Promotes Cell Apoptosis by Inducing Pendrin Transfer, Reducing Cl- Transport, and Inhibiting PI3K/Akt/mTOR Pathway *Biomed Res Int*. 2022 Aug 29;2022:6496799.

Version 2024.03.12