



碧云天生物技术/Beyotime Biotechnology
订货热线：400-168-3301或800-8283301
订货e-mail：order@beyotime.com
技术咨询：info@beyotime.com
网址：http://www.beyotime.com

H-89 (PKA抑制剂)

产品编号	产品名称	包装
S1643-1mg	H-89 (PKA抑制剂)	10mg/ml×0.1ml
S1643-5mg	H-89 (PKA抑制剂)	5mg
S1643-25mg	H-89 (PKA抑制剂)	25mg
S1643-100mg	H-89 (PKA抑制剂)	100mg

产品简介：

- H-89也称H89，全称为N-[2-(p-Bromocinnamylamino)ethyl]-5-isoquinolinesulfonamide · 2HCl hydrate，是一种常用的PKA抑制剂。H-89可以通透细胞，可以选择性抑制cAMP、cGMP依赖的蛋白激酶PKA、PKG和PKC μ ，但对于PKA的选择性抑制作用更强。
- H-89分子量为519.28，分子式为C₂₀H₂₀BrN₃O₂S · 2HCl，CAS Number: 127243-85-0。本产品纯度大于98%。
- 本H-89为进口分装，其中10mg/ml包装产品用DMSO配制，浓度接近20mM，共0.1ml。5mg, 25mg和100mg包装为粉末装。

包装清单：

产品编号	产品名称	包装
S1643-1mg	H-89 (PKA抑制剂)	10mg/ml×0.1ml
S1643-5mg	H-89 (PKA抑制剂)	5mg
S1643-25mg	H-89 (PKA抑制剂)	25mg
S1643-100mg	H-89 (PKA抑制剂)	100mg
—	说明书	1份

保存条件：

-20°C保存，一年有效。

注意事项：

- 本产品对人体有害，操作时请小心，并注意有效防护以避免直接接触人体或吸入体内。
- 本H-89在4°C、冰浴等较低温度情况下会凝固而粘在离心管管底、管壁或管盖内，可以20-25°C水浴温育片刻至全部融解后使用。
- 本产品仅限于专业人员的科学研究用，不得用于临床诊断或治疗，不得用于食品或药品，不得存放于普通住宅内。
- 为了您的安全和健康，请穿实验服并戴一次性手套操作。

使用说明：

1. H-89常见使用浓度范围为10-30 μ M，也有报道在1-50 μ M范围内使用。具体的最佳工作浓度请参考相关文献，或根据实验目的，以及所培养的特定细胞和组织，通过实验进行摸索和优化。

使用本产品的文献：

1. Yu Y, Liu L, Wang X, Liu X, Liu X, Xie L, Wang G. Modulation of glucagon-like peptide-1 release by berberine: in vivo and in vitro studies. Biochem Pharmacol. 2010;79(7):1000-6.
2. Li Y, Kang Z, Li S, Kong T, Liu X, Sun C. Ursolic acid stimulates lipolysis in primary-cultured rat adipocytes. Mol Nutr Food Res. 2010 Nov;54(11):1609-17.
3. Jiang Z, Li S, Liu Y, Deng P, Huang J, He G. Sesamin induces melanogenesis by microphthalmia-associated transcription factor and tyrosinase-up-regulation via cAMP signaling pathway. Acta Biochim Biophys Sin (Shanghai). 2011 Oct;43(10):763-70.
4. Zhang Q, Lu YQ, Jiang JK, Gu LH, Mou HZ. Early changes of CD4+CD25+Foxp3+ regulatory T cells and Th1/Th2, Tc1/Tc2 profiles in the peripheral blood of rats with controlled hemorrhagic shock and no fluid resuscitation. Chin Med J (Engl). 2012 Jun;125(12):2163-7.
5. Li J, Li Z, Tang L, Yang Y, Zouine M, Bouzayen M. A conserved phosphorylation site regulates the transcriptional function of ETHYLENE-INSENSITIVE3-like1 in tomato. J Exp Bot. 2012 Jan;63(1):427-39.
6. Liu Y, Su Y, Sun S, Wang T, Qiao X, Li H, Run X, Liang Z. Human tau may modify glucocorticoids-mediated regulation of cAMP-dependent kinase and phosphorylated cAMP response element binding protein. Neurochem Res. 2012 May;37(5):935-47.
7. Li Y, He J, Sui S, Hu X, Zhao Y, Li N. Clenbuterol upregulates histone demethylase JHDM2a via the β (2)-adrenoceptor/cAMP/PKA/p-CREB signaling pathway. Cell Signal. 2012 Dec;24(12):2297-306.
8. Mu X, He K, Sun H, Zhou X, Chang L, Li X, Chu W, Qiao G, Lu Y. Hydrogen peroxide induces overexpression of angiotensin-converting enzyme in human umbilical vein endothelial cells. Free Radic Res. 2013 Feb;47(2):116-22.
9. Chen Y, Su Y, Run X, Sun Z, Wang T, Sun S, Liang Z. Pretreatment of PC12 Cells with 17 β -estradiol Prevents A β -Induced Down-Regulation of CREB Phosphorylation and Prolongs Inhibition of GSK-3 β . J Mol Neurosci. 2013 Jul;50(3):394-401.
10. Chen ML, Yi L, Jin X, Liang XY, Zhou Y, Zhang T, Xie Q, Zhou X, Chang H, Fu YJ, Zhu JD, Zhang QY, Mi MT. Resveratrol attenuates vascular endothelial inflammation by inducing autophagy through the

cAMP signaling pathway. *Autophagy*. 2013 Dec;9(12):2033-45.

11. Ouyang C, Nie L, Gu M, Wu A, Han X, Wang X, Shao J, Xia Z. Transforming growth factor (TGF)- β -activated kinase 1 (TAK1) activation requires phosphorylation of serine 412 by protein kinase A catalytic subunit α (PKAC α) and X-linked protein kinase (PRKX). *J Biol Chem*. 2014 Aug 29;289(35):24226-37.
12. Du XH, Zhou XL, Cao R, Xiao P, Teng Y, Ning CB, Liu HL. FSH-induced p38-MAPK-mediated dephosphorylation at serine 727 of the signal transducer and activator of transcription 1 decreases Cyp1b1 expression in mouse granulosa cells. *Cell Signal*. 2015 Jan;27(1):6-14.
13. Wang Z, Zhang X, Tian H, Wang W, Ru S. Effects of monocrotophos pesticide on steroidogenesis and transcription of steroidogenic enzymes in rainbow trout RTG-2 cells involving the protein kinase A signal pathway. *Toxicol In Vitro*. 2015 Feb;29(1):155-61.
14. Zhang Y, Chen ML, Zhou Y, Yi L, Gao YX, Ran L, Chen SH, Zhang T, Zhou X, Zou D, Wu B, Wu Y, Chang H, Zhu JD, Zhang QY, Mi MT. Resveratrol improves hepatic steatosis by inducing autophagy through the cAMP signaling pathway. *Mol Nutr Food Res*. 2015 Aug;59(8):1443-57.
15. Wang J, Peng Y, Liu Y, Yang J, Huang M, Tan W. AT-101 inhibits hedgehog pathway activity and cancer growth. *Cancer Chemother Pharmacol*. 2015 Sep;76(3):461-9.
16. Chen C, Gu S, Jiang X, Zhang Z. Nuclear translocation of nuclear factor kappa B is regulated by G protein signaling pathway in arsenite-induced apoptosis in HBE cell line. *Environ Toxicol*. 2016 31: 1819-33.
17. Chen B, Lin T, Yang X, Li Y, Xie D, Cui H. Intermittent parathyroid hormone (1-34) application regulates cAMP-response element binding proteinactivity to promote the proliferation and osteogenic differentiation of bone mesenchymal stromal cells, via the cAMP/PKA signaling pathway. *Exp Ther Med*. 2016 Jun;11(6):2399-2406.
18. Shen C, Li X, Rasooly A, Guo L, Zhang K, Yang M. A single electrochemical biosensor for detecting the activity and inhibition of both protein kinase and alkaline phosphatase based on phosphate ions induced deposition of redox precipitates. *Biosens Bioelectron*. 2016 Nov 15;85:220-5.
19. Yu ZL, Li DQ, Huang XY, Xing X, Yu RQ, Li Z, Li ZB. Lysophosphatidic acid upregulates connective tissue growth factor expression in osteoblasts through the GPCR/PKC and PKA pathways. *Int J Mol Med*. 2016 Feb;37(2):468-74.
20. Luo QQ, Qian ZM, Zhou YF, Zhang MW, Wang D, Zhu L, Ke Y. Expression of Iron Regulatory Protein 1 Is Regulated not only by HIF-1 but also pCREB under Hypoxia. *Int J Biol Sci*. 2016 Sep 15;12(10):1191-1202.
21. Guo X, Yu R, Xu Y, Lian R, Yu Y, Cui Z, Ji Q, Chen J, Li Z, Liu H, Chen J. PAC1R agonist maxadilan enhances hADSC viability and neural differentiation potential. *J Cell Mol Med*. 2016 May;20(5):874-90.
22. Zhao J, Ma J, Lu J, Jiang Y, Zhang X, Zhao J, Yang H, Huang Y, Zhao M, Liu K, Dong Z. Involvement of p38MAPK-ATF2 signaling pathway in alternariol induced DNA polymerase β expression. *Oncol Lett*. 2016 Jul;12(1):675-679.
23. Meng M, Liao H, Zhang B, Pan Y, Kong Y, Liu W, Yang P, Huo Z, Cao Z, Zhou Q. Cigarette smoke extracts induce overexpression of the proto-oncogenic gene interleukin-13 receptor $\alpha 2$ through activation of the PKA-CREB signaling pathway to trigger malignant transformation of lung vascular endothelial cells and angiogenesis. *Cell Signal*. 2017 Feb;31:15-25.

Version 2016.12.09