



碧云天生物技术/Beyotime Biotechnology
订货热线：400-168-3301或800-8283301
订货e-mail：order@beyotime.com
技术咨询：info@beyotime.com
网址：http://www.beyotime.com

MG-132 (Proteasome抑制剂)

产品编号	产品名称	包装
S1748-1mg	MG-132 (Proteasome抑制剂)	20mg/ml×0.05ml
S1748-5mg	MG-132 (Proteasome抑制剂)	5mg
S1748-25mg	MG-132 (Proteasome抑制剂)	25mg

产品简介：

- MG-132也写作MG132，也称Z-LLL-CHO、Z-Leu-Leu-Leu-CHO、Z-Leu-Leu-Leu-al或Carbobenzoxy-L-leucyl-L-leucyl-L-leucinal，是一种常用的Proteasome抑制剂，即蛋白酶体抑制剂。MG-132可以通透细胞，选择性抑制Proteasome，Ki=4nM。可以激活JNK1并抑制NF-κB的激活(IC50=3μM)。在10μM时可以诱导PC12细胞产生neurite outgrowth。
- MG-132分子量为475.62，分子式为C₂₆H₄₁N₃O₅，CAS Number: 133407-82-6。本产品纯度大于99%。
- 本MG-132为进口分装，其中20mg/ml包装产品用DMSO配制，共0.05ml。5mg和25mg包装为粉末装。

包装清单：

产品编号	产品名称	包装
S1748-1mg	MG-132 (Proteasome抑制剂)	20mg/ml×0.05ml
S1748-5mg	MG-132 (Proteasome抑制剂)	5mg
S1748-25mg	MG-132 (Proteasome抑制剂)	25mg
—	说明书	1份

保存条件：

-20°C避光保存，半年有效。

注意事项：

- 本MG-132在4°C、冰浴等较低温度情况下会凝固而粘在离心管管底、管壁或管盖内，可以20-25°C水浴温育片刻至全部融解后使用。
- 本产品仅限于专业人员的科学研究用，不得用于临床诊断或治疗，不得用于食品或药品，不得存放于普通住宅内。
- 为了您的安全和健康，请穿实验服并戴一次性手套操作。

使用说明：

1. MG-132常见使用浓度范围为1-50μM。具体的最佳工作浓度请参考相关文献，或根据实验目的，以及所培养的特定细胞和组织，通过实验进行摸索和优化。

使用本产品的文献：

1. Wu H, Chen Y, Wang ZY, Li W, Li JQ, Zhang L, Lu YJ. Involvement of p21 (waf1) in merlin deficient sporadic vestibular schwannomas. Neuroscience. 2010;170(1):149-55.
2. Wen L, Liu J, Chen Y, Wu D. Identification and preliminary functional analysis of alternative splicing of Siah1 in Xenopus laevis. Biochem Biophys Res Commun. 2010;396(2):419-24.
3. Li L, Cui D, Zheng SJ, Lou H, Tang J. Regulation of Actinomycin D induced upregulation of Mdm2 in H1299 cells. DNA Repair (Amst). 2012 Feb 1;11(2):112-9.
4. Huang HL, Weng HY, Wang LQ, Yu CH, Huang QJ, Zhao PP, Wen JZ, Zhou H, Qu LH. Triggering Fbw7-mediated proteasomal degradation of c-Myc by oridonin induces cell growthinhibition and apoptosis. Mol Cancer Ther. 2012 May;11(5):1155-65.
5. Wang D, Fang L, Li K, Zhong H, Fan J, Ouyang C, Zhang H, Duan E, Luo R, Zhang Z, Liu X, Chen H, Xiao S. Foot-and-Mouth Disease Virus 3C Protease Cleaves NEMO To Impair Innate Immune Signaling. J Virol. 2012 Sep;86(17):9311-22.
6. Lu L, Song HF, Zhang WG, Liu XQ, Zhu Q, Cheng XL, Yang GJ, Li A, Xiao ZC. Potential role of 20S proteasome in maintaining stem cell integrity of human bone marrow stromalcells in prolonged culture expansion. Biochem Biophys Res Commun. 2012 May 25;422(1):121-7.
7. Wang L, Chen M, Yang J, Zhang Z. LC3 fluorescent puncta in autophagosomes or in protein aggregates can be distinguished by FRAP analysis inliving cells. Autophagy. 2013 May;9(5):756-69.
8. Qiu M, Chen Y, Chu Y, Song S, Yang N, Gao J, Wu Z. Zinc ionophores pyritohone inhibits herpes simplex virus replication through interfering with proteasome functionand NF-κB activation. Antiviral Res. 2013 Oct;100(1):44-53.
9. Chen HY, Ren XY, Wang WH, Zhang YX, Chen SF, Zhang B, Wang LX. Upregulated ROS production induced by the proteasome inhibitor MG-132 on XBPI gene expression and cell apoptosis in Tca-8113 cells. Biomed Pharmacother. 2014 Jul;68(6):709-13.
10. Zhou Y, Chen S, Ding R, Yao W, Gao X. Inflammatory modulation effect of glycopeptide from Ganoderma capense (Lloyd) Teng. Mediators Inflamm. 2014;2014:691285.

11. Liu Z, Chen X, Wang Y, Peng H, Wang Y, Jing Y, Zhang H. PDK4 protein promotes tumorigenesis through activation of cAMP-response element-binding protein (CREB)-Ras homolog enriched in brain (RHEB)-mTORC1 signaling cascade. *J Biol Chem.* 2014 Oct 24;289(43):29739-49.
12. Liang Y, Hu J, Li J, Liu Y, Yu J, Zhuang X, Mu L, Kong X, Hong D, Yang Q, Hu G. Epigenetic Activation of TWIST1 by MTDH Promotes Cancer Stem-like Cell Traits in Breast Cancer. *Cancer Res.* 2015 Sep 1;75(17):3672-80.
13. Dong QM, Ling C, Zhu JF, Chen X, Tang Y, Zhao LI. Correlation between p65 and TNF- α in patients with acute myelocytic leukemia. *Oncol Lett.* 2015 Nov;10(5):3305-3309. 2015 Nov;10(5):3305-3309.
14. Xu Y, Liu F, Liu J, Wang D, Yan Y, Ji S, Zan J, Zhou J. The co-chaperone Cdc37 regulates the rabies virus phosphoprotein stability by targeting to Hsp90AA1 machinery. *Sci Rep.* 2016 Jun 2;6:27123.
15. Zhang C, Yang L, Geng YD, An FL, Xia YZ, Guo C, Luo JG, Zhang LY, Guo QL, Kong LY. Icariside II, a natural mTOR inhibitor, disrupts aberrant energy homeostasis via suppressing mTORC1-4E-BP1 axis in sarcoma cells. *Oncotarget.* 2016 May 10;7(19):27819-37.
16. Chen X, Cao X, Sun X, Lei R, Chen P, Zhao Y, Jiang Y, Yin J, Chen R, Ye D, Wang Q, Liu Z, Liu S, Cheng C, Mao J, Hou Y, Wang M, Siebenlist U, Eugene Chin Y, Wang Y, Cao L, Hu G, Zhang X. Bcl-3 regulates TGF β signaling by stabilizing Smad3 during breast cancer pulmonary metastasis. *Cell Death Dis.* 2016 Dec 1;7(12):e2508.

Version 2016.12.12